

# Sterilized filter bag for Homogenize PYXON-20 *sereise*

(Worldwide PATENT Pending)

&lt;MORE THAN LESS LEES&gt; &lt;EASY TO USE PIPET&gt; &lt;DISPENSER PIPETTE USABLE&gt; &lt;INEXPENSIVE&gt;

## WIDE MOUTH

Easier to open with a wide filter opening ! Sample inlet is also wide.

## DISPENSER (Micro pipette) CAN BE USED

### MAKE A SLIT

For dispenser (Micro pipette) use, slit the bag with picking Pull-Tab and down until Pull down guide line where is printed at the left bottom for pull down aim.

#### Pull down guide lines

The upper line is for 250ml

The lower line is for 100ml

In case of 10g specimen and 10 times sterilized water added, the water level will be about 4cm under the lower line.

#### Be careful

Do not slit down under the line. Slit straight when you slit pull-tab down.

### DISPENSER OPERATION

Insert your Dispenser from the Slit made as the above way, then, suck specimen up.



## WRITE-ON

It's easy to write data on the white print with an ordinary ballpoint pen, pencil, or felt-tip marker.

## SANITARY PORTION PACK

All items are packed by zip bag.

### SPECIFICATIONS

Product name	Code No.	Specifications	Qty/box	Bag size	Capacity
Sterilized homogenize bag	PX0010	Plain homogenize bag without Filter.	1,000pcs	180×300mm	400ml
Sterilized filter bag	PX0020	Pull-down type for a micro-pipette use	500pcs	190×300mm	400ml
	PX0020P	Plain type for a measuring pipette use	500pcs	190×300mm	400ml
	PX0020S	For Spiral plating method use	500pcs	190×300mm	400ml
	PX0020M	Small size filter bag with pull-down tab	1,000pcs	114×190mm	80ml

Specifications are subject to change without notice for remodeling or quality enhancement.

Manufacturer

**ELMEX Limited**

2-2, Ichigayasadohara-cho, Shinjuku-ku, Tokyo 162-0842 Japan

Tel: 81-3-5229-7911 Fax: 81-3-5261-4130

URL://www.elmex.co.jp e-mail: sales@elmex.co.jp